DPA-LOP.002 Attachment C

| Date | Analyst | | | | |
|-------------------------------|---------|---------|----------|-----------------------------------------|--|
| Dissolution Apparatus: Manufa | acturer | Model # | Serial # | Glass Dissolution Vessels: Manufacturer | |

MECHANICAL CALIBRATION REPORT SHEET -- APPARATUS 2 (PADDLE)

| Calibration | Point of | Results & Comments | Tools | DPA |
|---------------------------------------------------------------------------------|------------------------------------------------------------------------------------------------------------------------|---------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|----------------------------|---------------------------------------------------------------------------------------------|
| Parameter | Measurement | | Used | Tolerances |
| Vessel dimensions (Inspect at time of receipt only) | See USP | | Caliper, ruler | See USP |
| Paddle dimensions (Inspect at time of receipt only) | See USP | | Caliper, ruler | See USP |
| Paddle examination (Examine at time of use) | | | | No gross defects as determined by visual inspection |
| Visual inspection of belts, check ball bearings, oil bearings and rollers | | | | Belts are tight, clean, free of cracks, properly aligned, shafts turn freely |
| Shaft wobble | 2 cm above top of paddle blade | 1 2 3 4 5 6 | Runout gauge | \leq 0.5 mm total runout |
| Shaft verticality | Above vessel | Record results at 2 points that are 90° apart. Bubble is centered within lines at the following positions: (Y/N) Shaft1 Pt1: Pt2: Shaft4 Pt1: Pt2: Shaft2 Pt1: Pt2: Shaft5 Pt1: Pt2: Shaft3 Pt1: Pt2: Shaft6 | Bubble level | Bubble centered within the lines of a bubble level |
| Vessel/Shaft centering and vessel verticality | Setp 1: A Distek CenterChek 1.5 mm above blade. Step 2: Move CenterChek 80 mm above blade. Note: Blade bottom 2.5 cm | Step 1: CenterChek 1.5 mm above blade 1. 2. 3. 4. 5. 6. Step 2: CenterChek 80 mm above blade 1. 2. 3. 4. 5. 6. | Distek CenterChek | 1 mm from centerline |
| Height check/Paddle depth | above vessel bottom Paddle Bottom | 1 2 3 4 5 6 | Distek HeightChek | 25 ± 2 mm |
| Vibration | Paddle shafts, baseplete | | None | No significant vibration (tolerance may be changed upon further vibration studies) |
| Rotational speed | | 50 rpm 100 rpm | Digital optical tachometer | ± 2 rpm |
| Vessel temperature (Measured at time of use) | | 1 2 3 4 5 6 | Mercury thermometer | 37 ± 0.5° C |